# EASYCAL™ 5

Server Installation

# Impressum

BRAND GMBH + CO KG  
Otto-Schott-Str. 25  
97877 Wertheim (Germany)  
T +49 9342 808 0  
F +49 9342 808 98000  
[info@brand.de](mailto:info@brand.de)  
[www.brand.de](http://www.brand.de)

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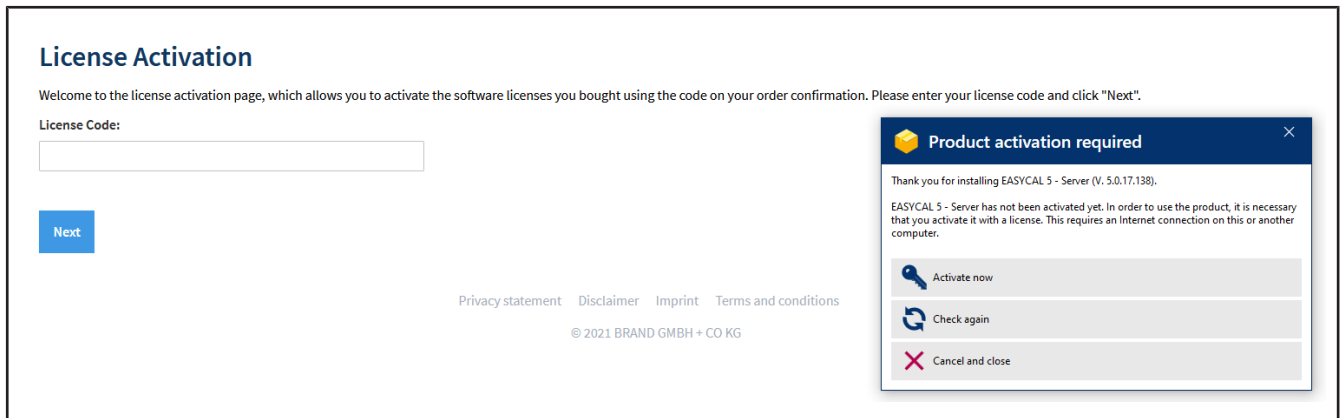
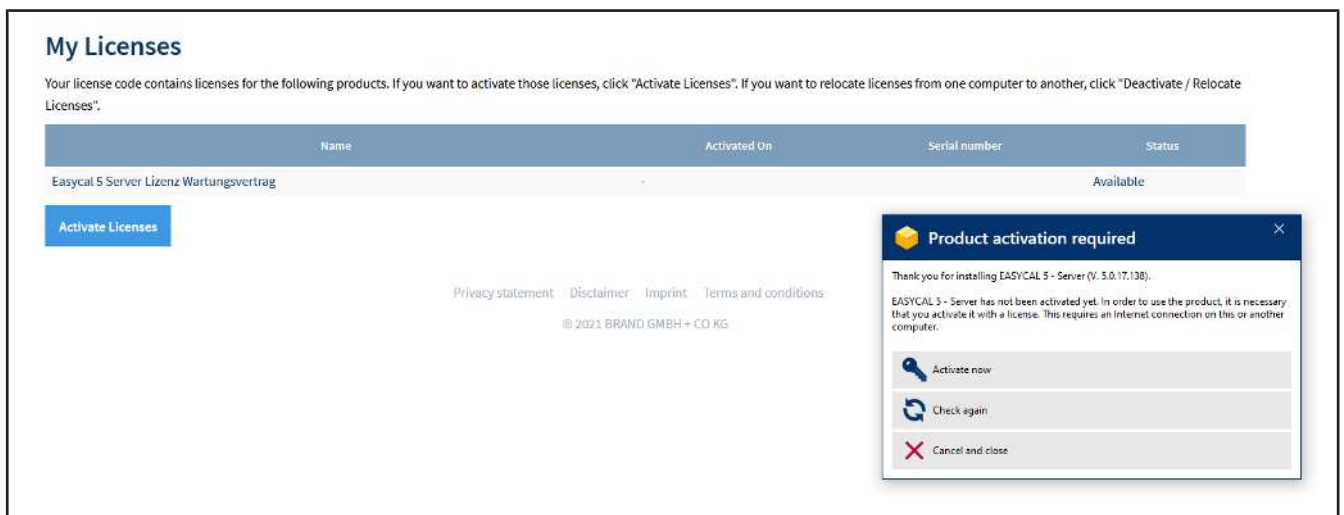
The screenshot shows the 'License Activation' page. At the top, it says 'License Activation' and 'Welcome to the license activation page, which allows you to activate the software licenses you bought using the code on your order confirmation. Please enter your license code and click "Next".' Below this is a 'License Code:' label and an empty text input field. A blue 'Next' button is positioned below the input field. At the bottom of the page, there are links for 'Privacy statement', 'Disclaimer', 'Imprint', and 'Terms and conditions', along with the copyright notice '© 2021 BRAND GMBH + CO KG'. A modal dialog box titled 'Product activation required' is overlaid on the right side. The dialog contains the text: 'Thank you for installing EASYCAL 5 - Server (V. 5.0.17.138). EASYCAL 5 - Server has not been activated yet. In order to use the product, it is necessary that you activate it with a license. This requires an Internet connection on this or another computer.' Below the text are three buttons: 'Activate now' (with a key icon), 'Check again' (with a refresh icon), and 'Cancel and close' (with a close icon).

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The screenshot shows the 'My Licenses' page. At the top, it says 'My Licenses' and 'Your license code contains licenses for the following products. If you want to activate those licenses, click "Activate Licenses". If you want to relocate licenses from one computer to another, click "Deactivate / Relocate Licenses".' Below this is a table with the following columns: 'Name', 'Activated On', 'Serial number', and 'Status'. The table contains one row with the following data: 'Easycal 5 Server Lizenz Wartungsvertrag', '-', and 'Available'. Below the table is a blue 'Activate Licenses' button. At the bottom of the page, there are links for 'Privacy statement', 'Disclaimer', 'Imprint', and 'Terms and conditions', along with the copyright notice '© 2021 BRAND GMBH + CO KG'. A modal dialog box titled 'Product activation required' is overlaid on the right side, identical to the one in the previous screenshot.

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[Activate now](#)  
[Check again](#)  
[Cancel and close](#)

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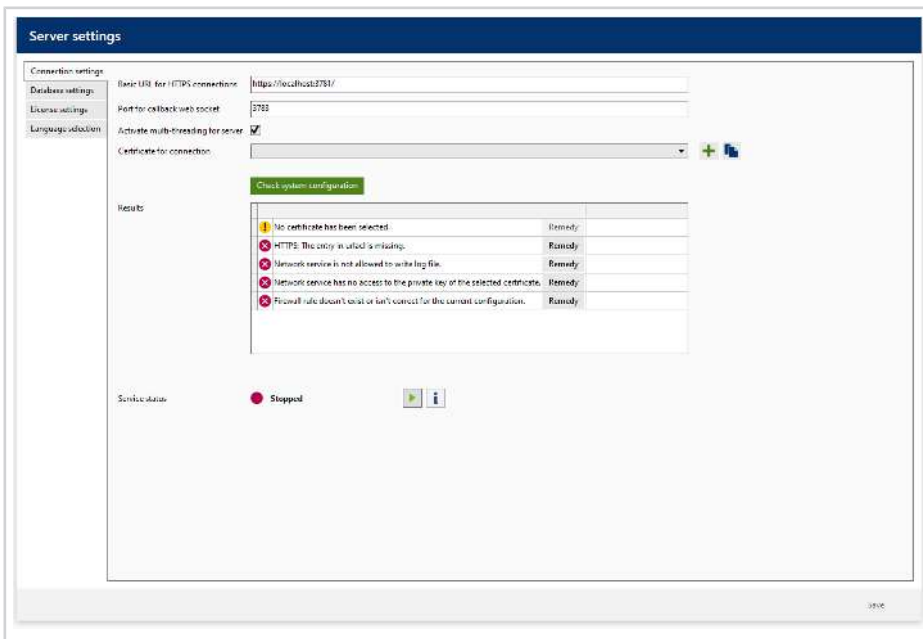
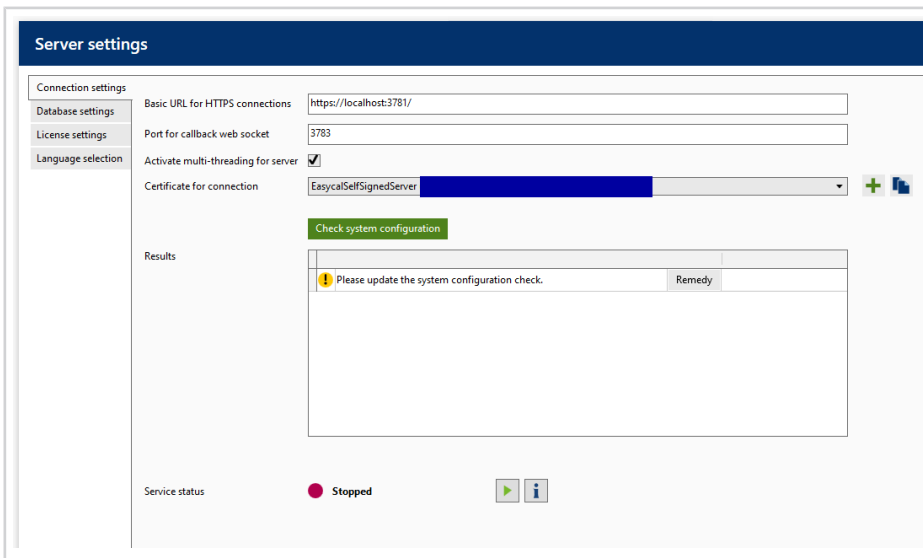
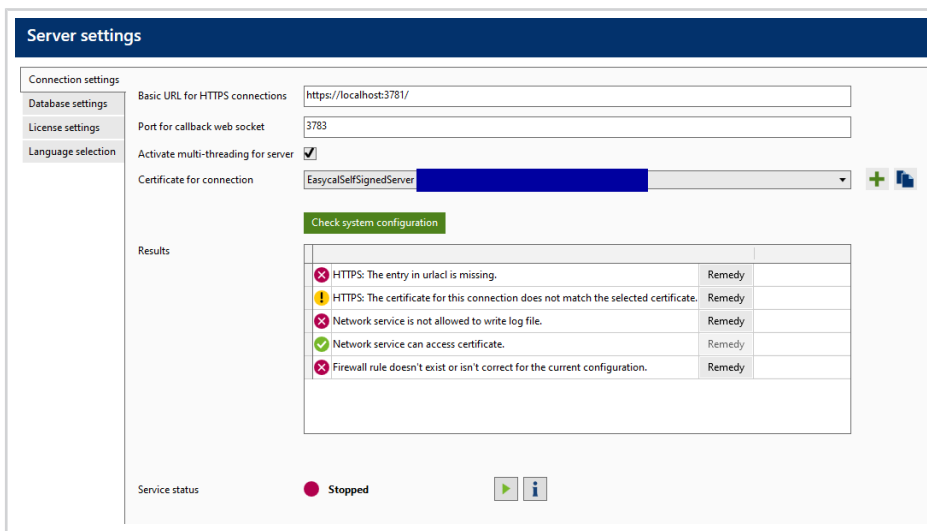
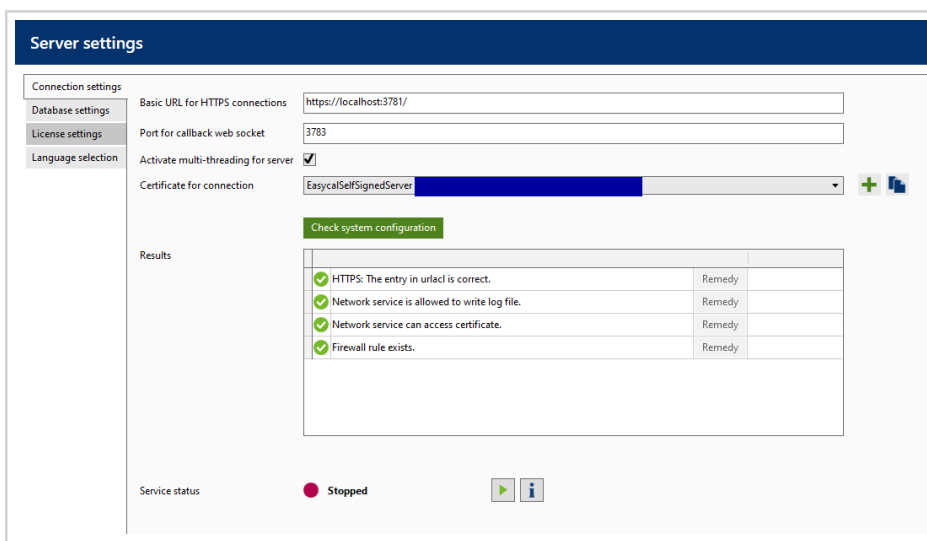
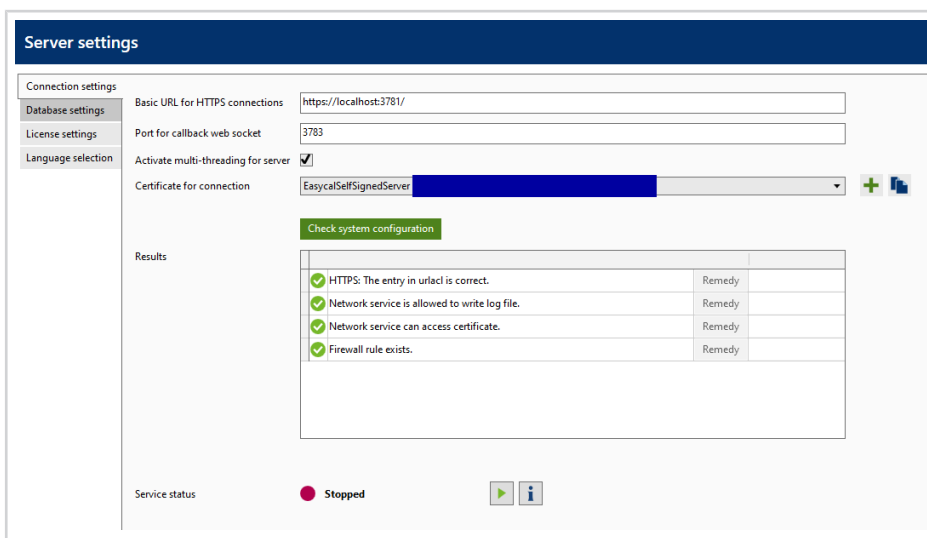
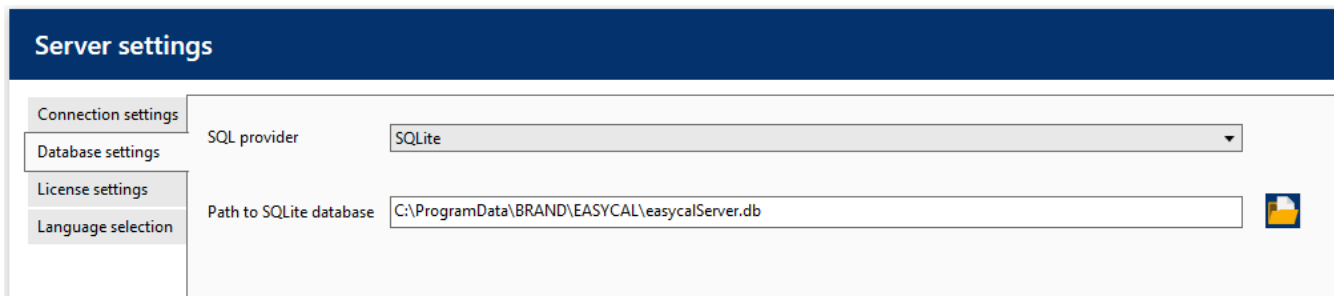
The screenshot shows the 'Server settings' dialog box with the 'Database settings' tab selected. The 'SQL provider' dropdown menu is set to 'SQLite'. The 'Path to SQLite database' text box contains the path 'C:\ProgramData\BRAND\EASYCAL\easycalServer.db'. There is a folder icon button to the right of the text box. On the left side, there are four tabs: 'Connection settings', 'Database settings', 'License settings', and 'Language selection'.

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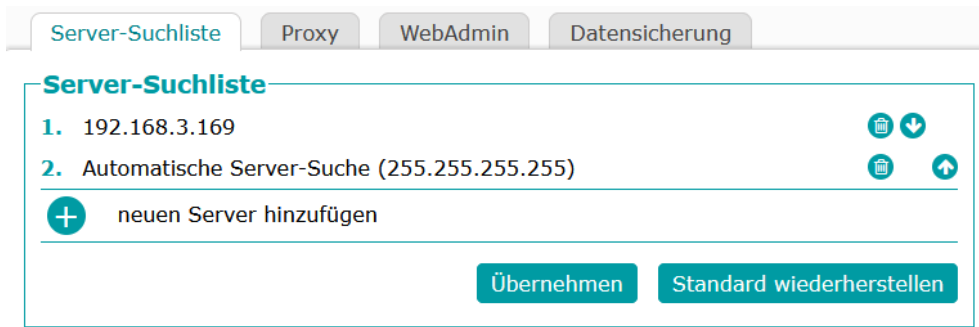
More operating manuals and translations?

Please refer to <http://www.brand.de/om> or use the following Quick Response Code:

Subject to technical changes, errors, and misprints.

# Contents

<b>1</b>	<b>System requirements.....</b>	<b>24</b>
<b>2</b>	<b>Installation .....</b>	<b>25</b>
<b>3</b>	<b>Activate License .....</b>	<b>27</b>
<b>4</b>	<b>Server software settings .....</b>	<b>30</b>
<b>5</b>	<b>Database settings .....</b>	<b>32</b>
<b>6</b>	<b>Add client licenses. ....</b>	<b>33</b>
<b>7</b>	<b>First Client Start .....</b>	<b>36</b>
<b>8</b>	<b>Configure the license server on the client .....</b>	<b>41</b>

# 1 System requirements

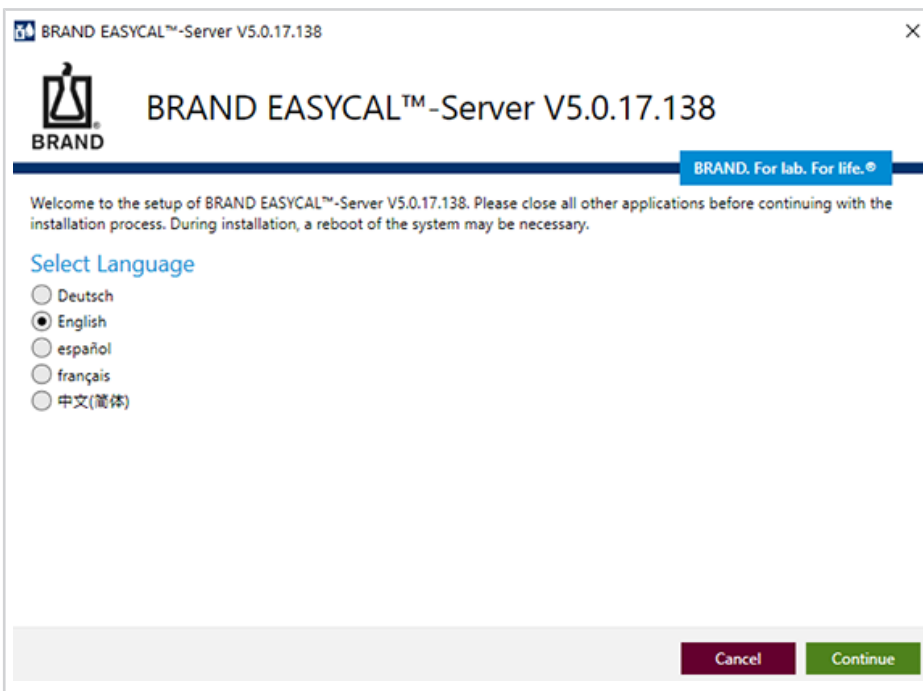
## System requirements for EASYCAL™ 5 as server

	<b>System requirements</b>
CPU:	Processor with at least 1 GHz clock rate
Memory	At least 16 GB RAM
Network	Local network infrastructure
Operating system	from Windows 10 or Windows Server 2016
Database	Local storage If a server operating system is used, the databases can also be stored locally on this operating system without having to use an SQL server. Alternatively: SQL Server as of MicrosoftSQL Server 2016
Internet	A permanent broadband Internet connection is highly recommended

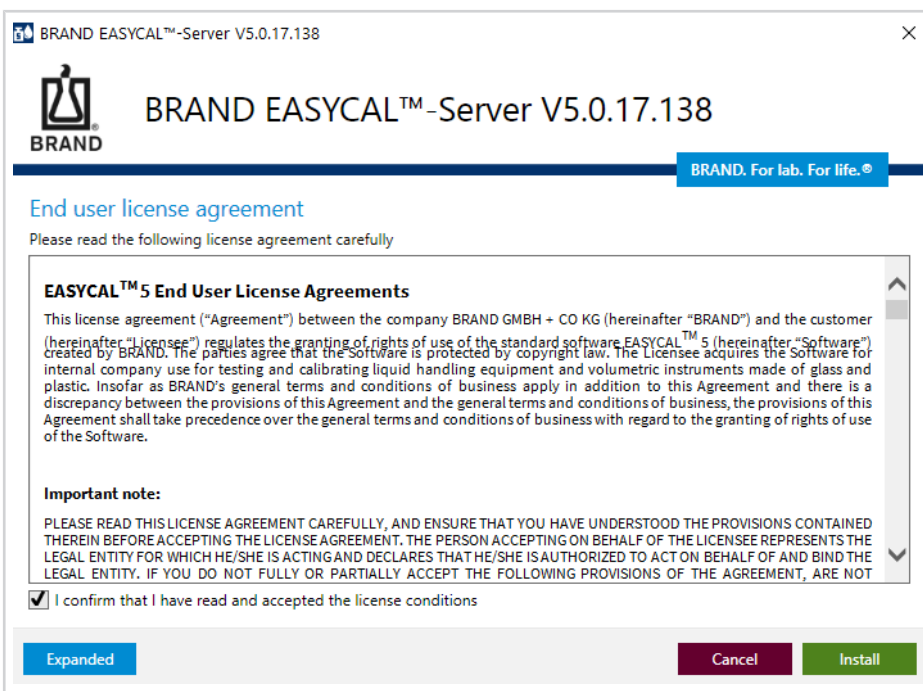
## 2 Installation

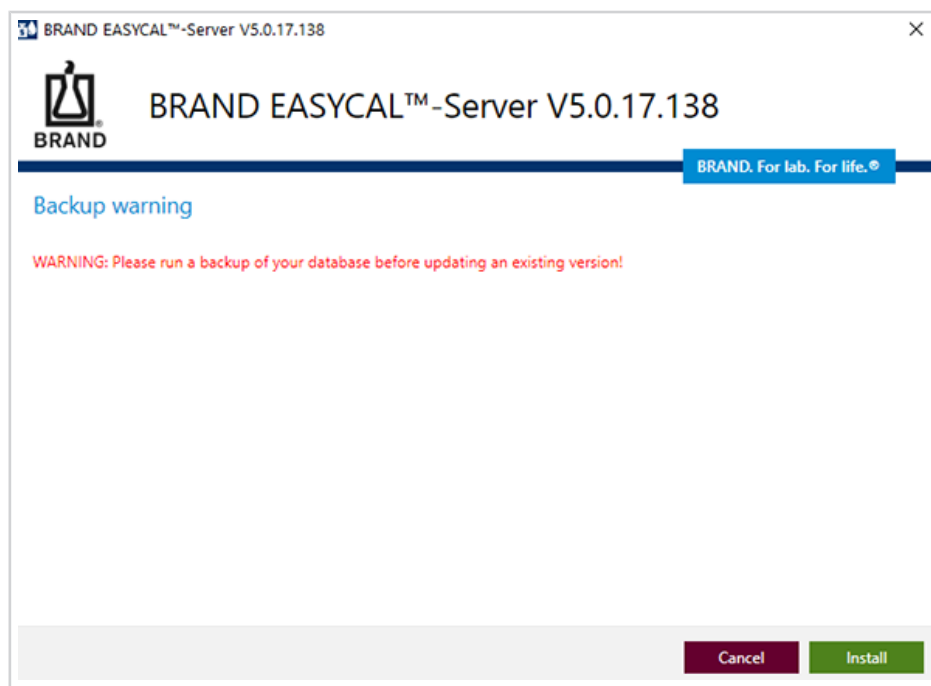
- a. Download the installation file: <https://update.brand.de/EASYCAL5Server/EASYCAL-Server-Setup-lat-est.exe>
- b. Start the installer using local admin rights and follow the dialog.
- c. Choose installation language. You can change the software language later on.

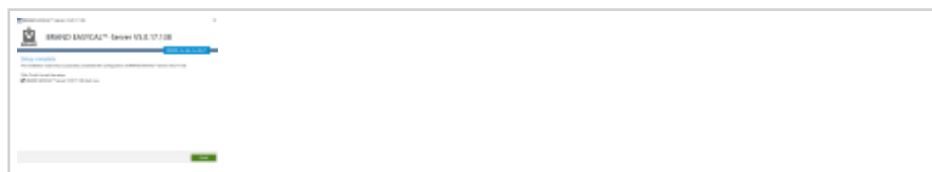
- d. Confirm the license agreement.

- e. If EASYCAL™ 5 has been installed before, we recommend backing up the database. Otherwise, you can proceed with the installation.

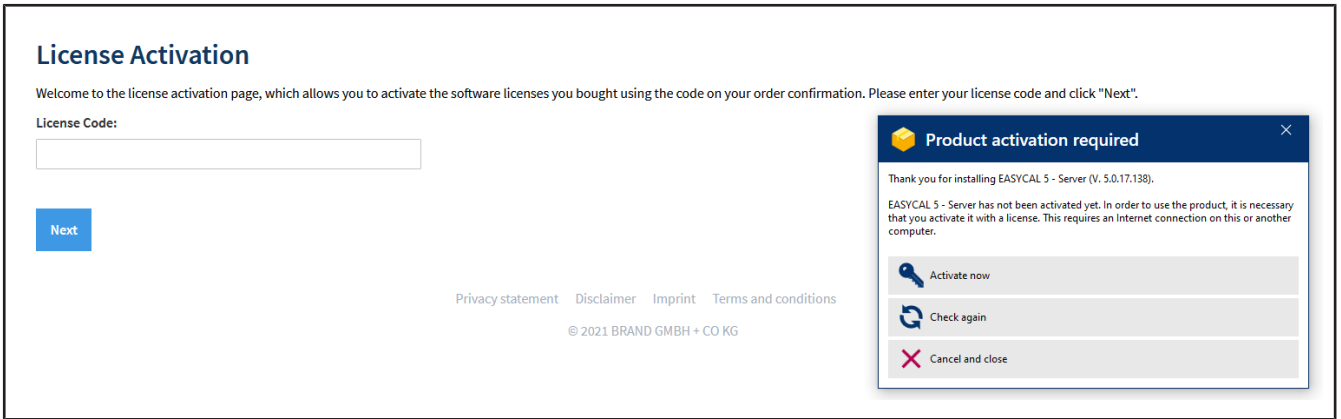
- f. Start the server after installation.

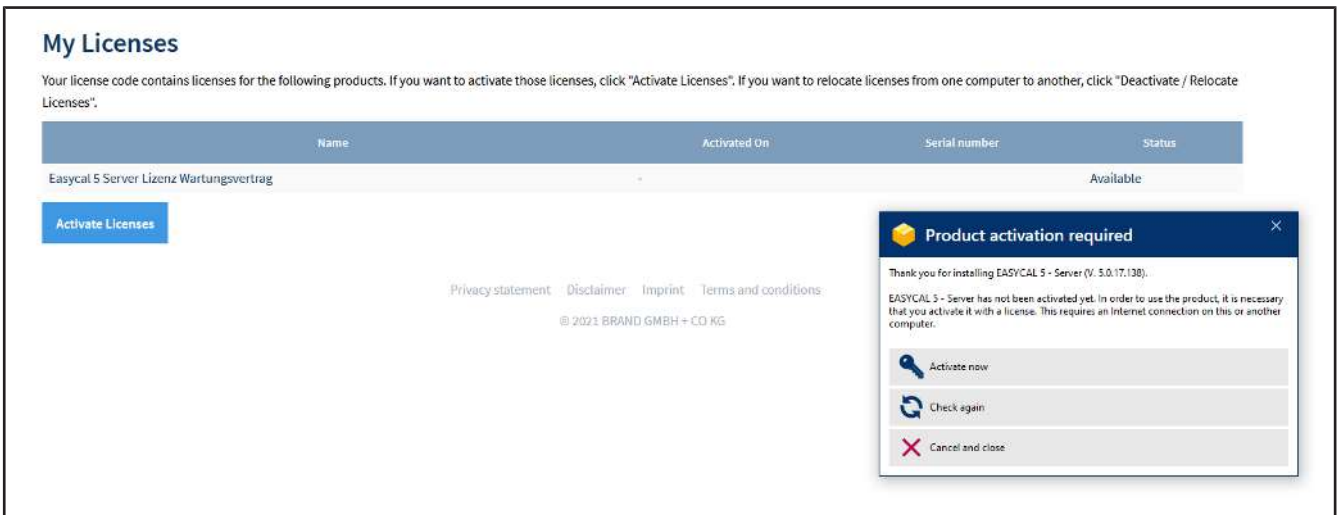
### 3 Activate License

When the EASYCAL™ 5 server software is started for the first time, the context menu for product activation opens.

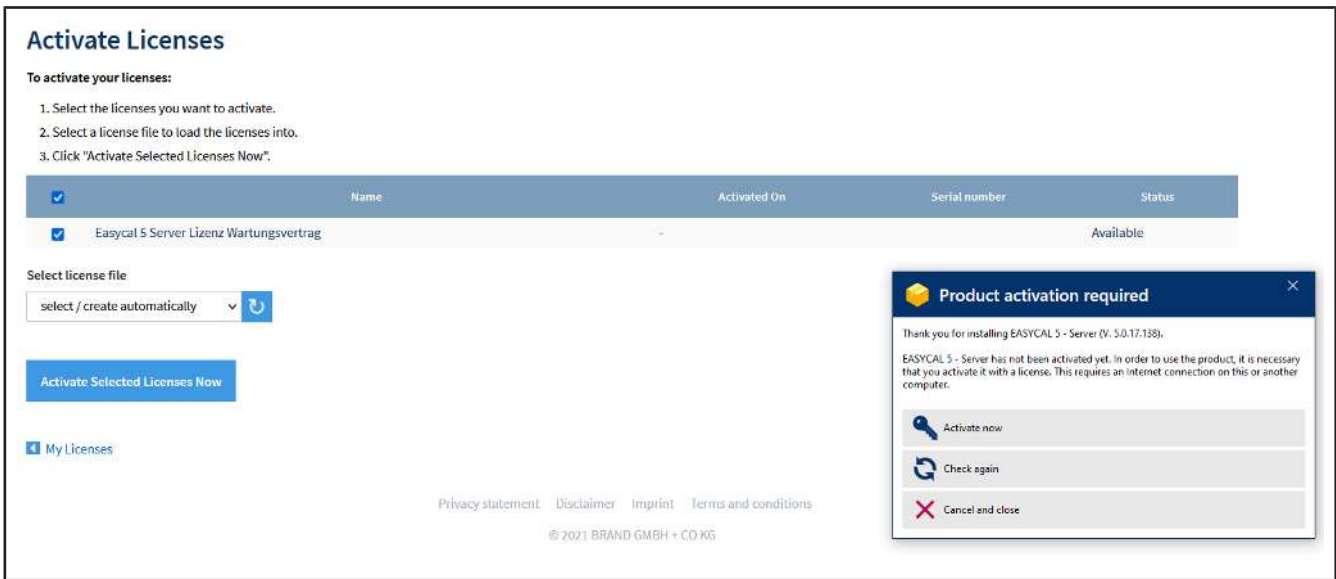
- a. Click Activate now, your default web browser will open.

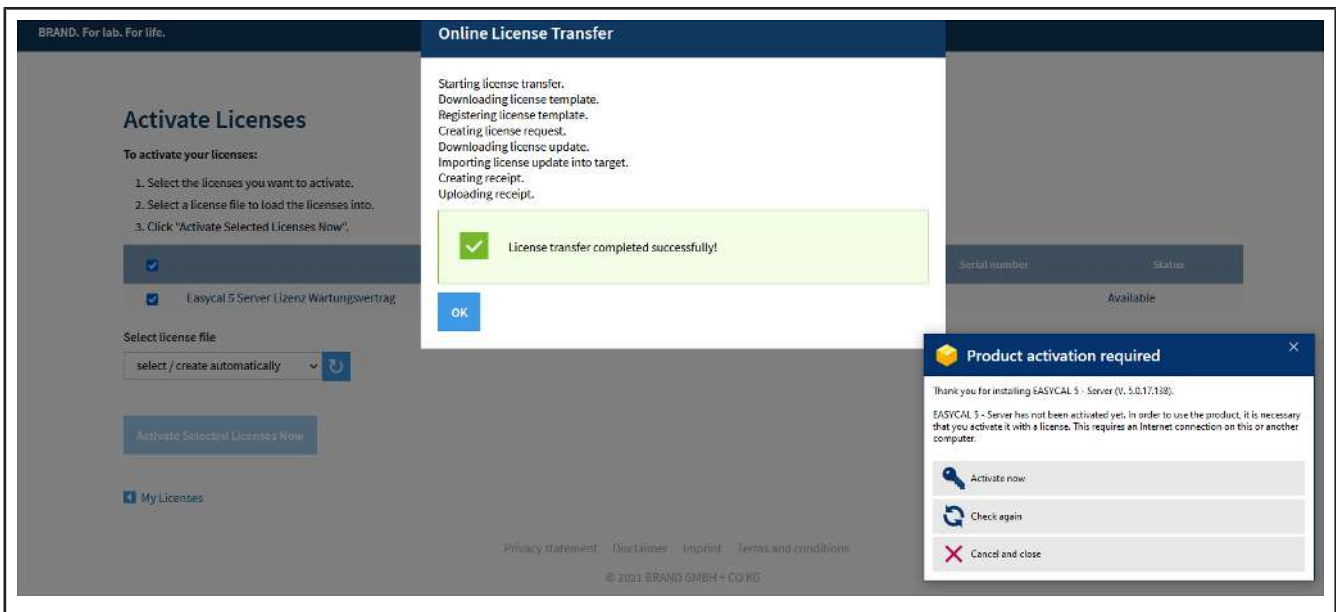
- b. Enter the license code for the EASYCAL™ 5 server software and click Next.

- c. Click on Activate Licenses.

- d. We recommend to keep the standard setting. Choose the EAYSCAL™ 5 Server License or the EAYSCAL™ 5 Server License Maintenance Contract - depending on which product you have purchased. Click on Activate Selected Licenses Now.

- e. The licensing process is performed. Confirm with OK.

**My Licenses**

Your license code contains licenses for the following products. If you want to activate those licenses, click "Activate Licenses". If you want to relocate licenses from one computer to another, click "Deactivate / Relocate Licenses".

Name	Activated On	Serial number	Status
Easycal 5 Server Lizenz Wartungsvertrag	2022-01-17 15:20:12	130-474324970	Activated

[Deactivate / Relocate Licenses](#)
[Restore Licenses](#)

[Privacy statement](#)
[Disclaimer](#)
[Imprint](#)
[Terms and conditions](#)

© 2021 BRAND GMBH + CO KG

**Product activation required** ✕

Thank you for installing EASYCAL 5 - Server (V. 5.0.17.138).

EASYCAL 5 - Server has not been activated yet. In order to use the product, it is necessary that you activate it with a license. This requires an Internet connection on this or another computer.

🔑 [Activate now](#)

🔄 [Check again](#)

✖ [Cancel and close](#)

f. In the Product Activation Required window, click Check Again.

**My Licenses**

Your license code contains licenses for the following products. If you want to activate those licenses, click "Activate Licenses". If you want to relocate licenses from one computer to another, click "Deactivate / Relocate Licenses".

Name	Activated On	Serial number	Status
Easycal 5 Server Lizenz Wartungsvertrag	2022-01-17 15:20:12	130-474324970	Activated

[Deactivate / Relocate Licenses](#)
[Restore Licenses](#)

[Privacy statement](#)
[Disclaimer](#)
[Imprint](#)
[Terms and conditions](#)

© 2021 BRAND GMBH + CO KG

**Product activation required** ✕

Thank you for installing EASYCAL 5 - Server (V. 5.0.17.138).

EASYCAL 5 - Server has not been activated yet. In order to use the product, it is necessary that you activate it with a license. This requires an Internet connection on this or another computer.

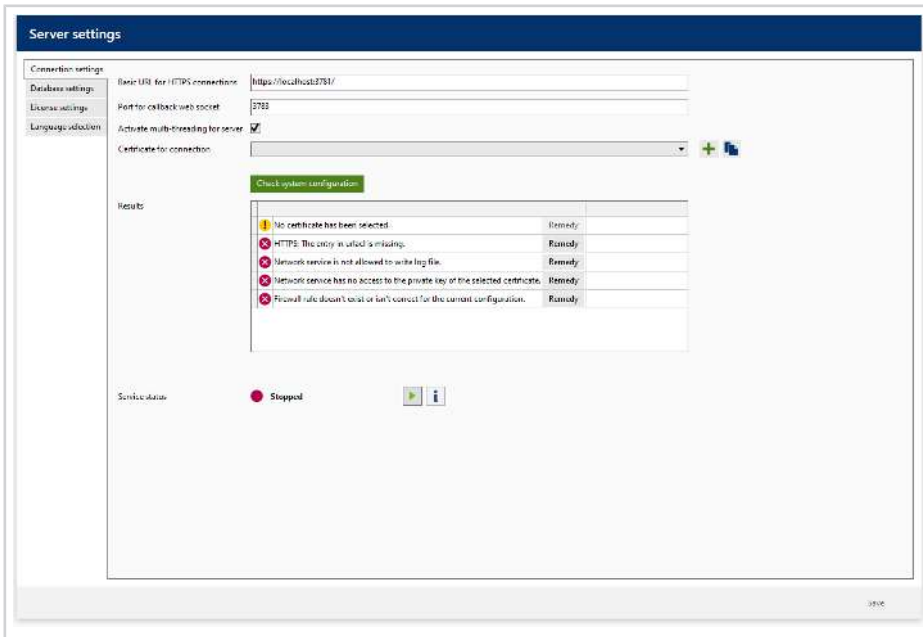
🔑 [Activate now](#)

🔄 [Check again](#)

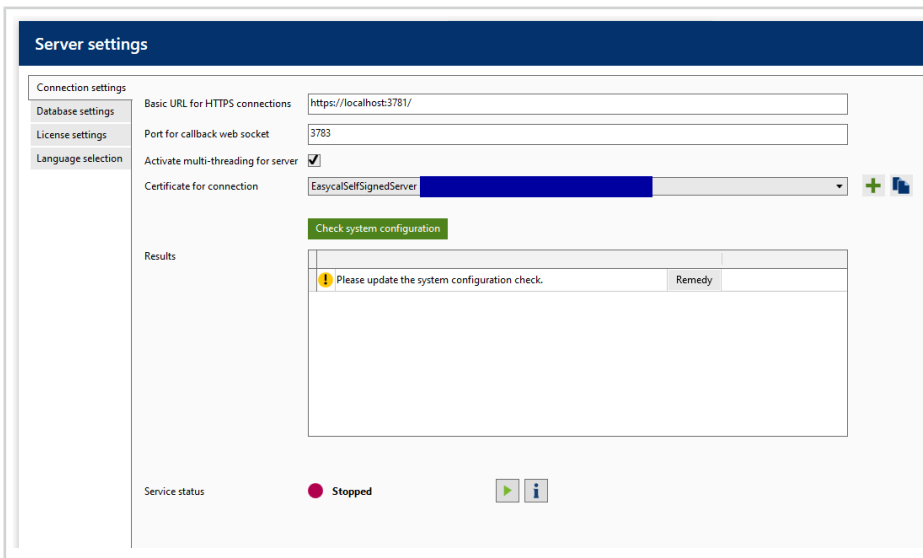
✖ [Cancel and close](#)

## 4 Server software settings

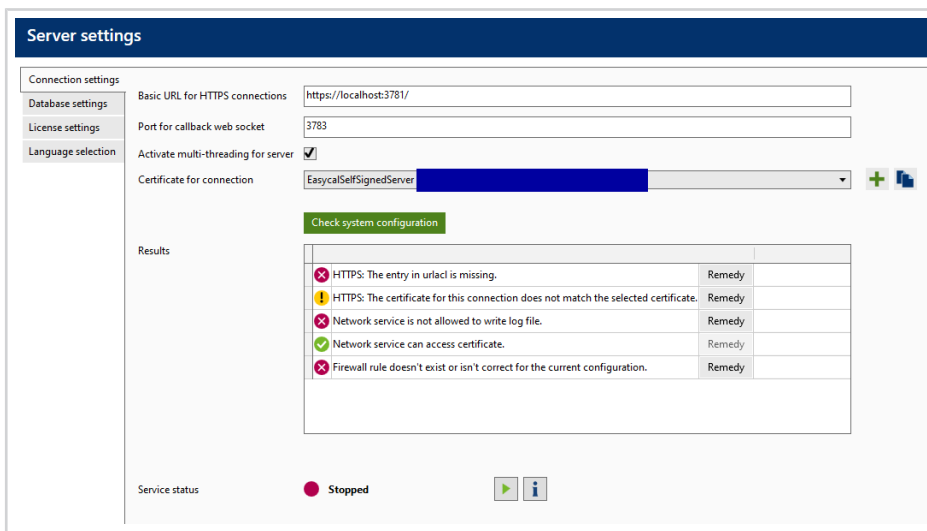
a. After clicking Check Again in the Product activation required window, the server software opens.

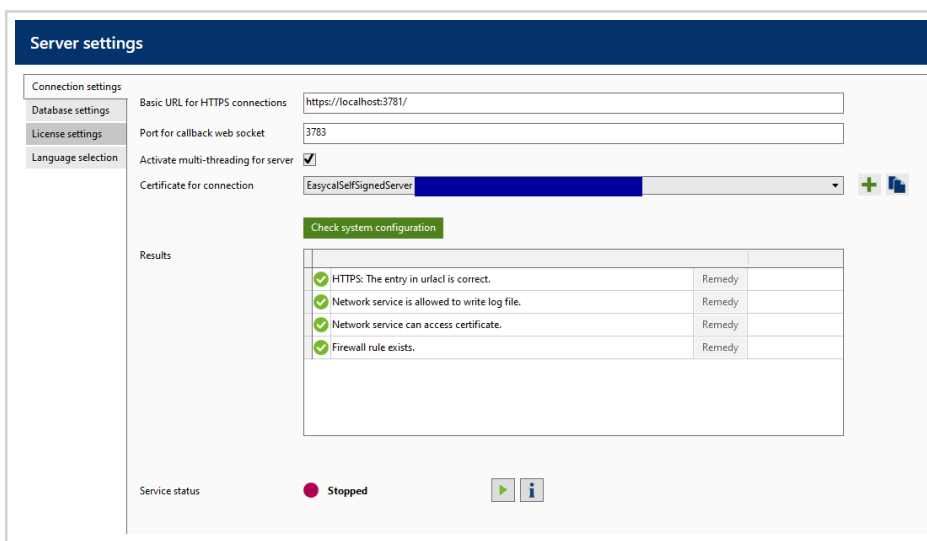
b. Click on the plus symbol at Certificate for connection and select EasycaSelfSignedServer.

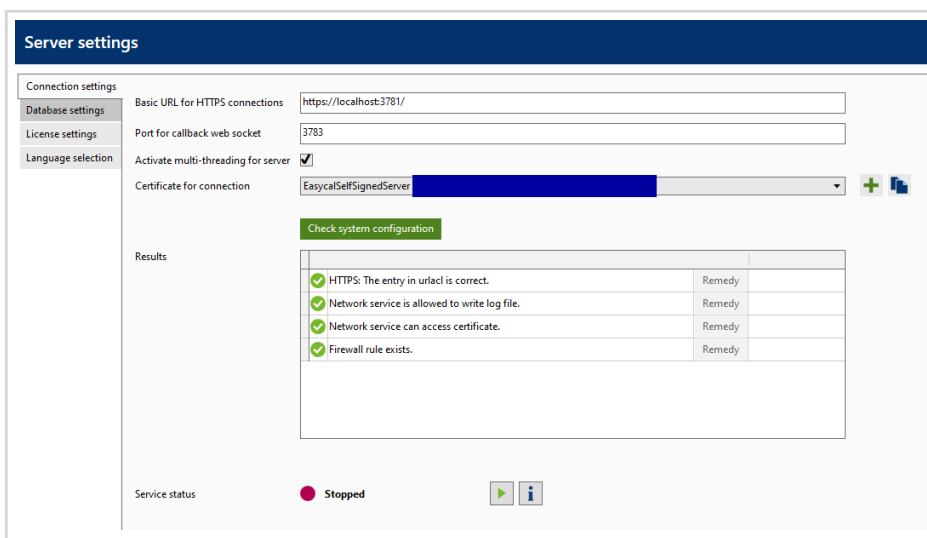
c. Then click on Check system configuration .

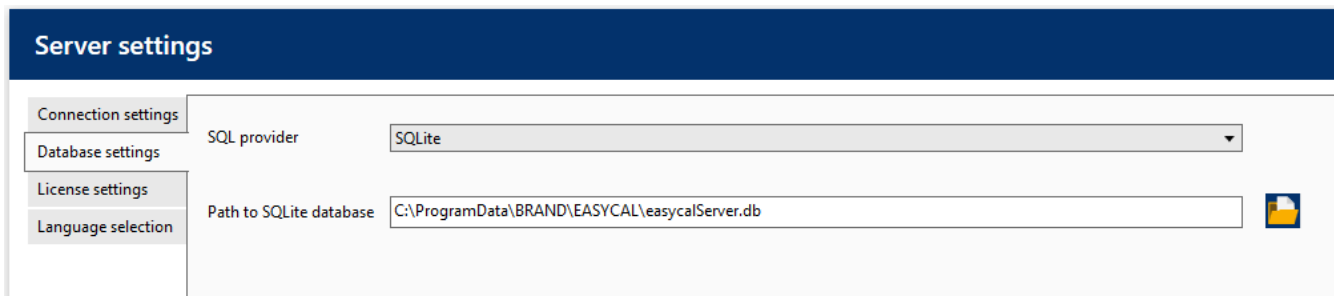
Click on Remedy for each of the Results so that the Cooling Block server software creates suitable entries in the system.

d. Click on Save.

## 5 Database settings

The screenshot shows a window titled "Server settings" with a dark blue header. On the left, there is a vertical sidebar with four tabs: "Connection settings", "Database settings", "License settings", and "Language selection". The "Database settings" tab is selected. The main area contains two settings:

- "SQL provider": A dropdown menu with "SQLite" selected.
- "Path to SQLite database": A text input field containing "C:\ProgramData\BRAND\EASYCAL\easyCalServer.db" and a folder icon button to the right.

Switch to the Database settings tab.

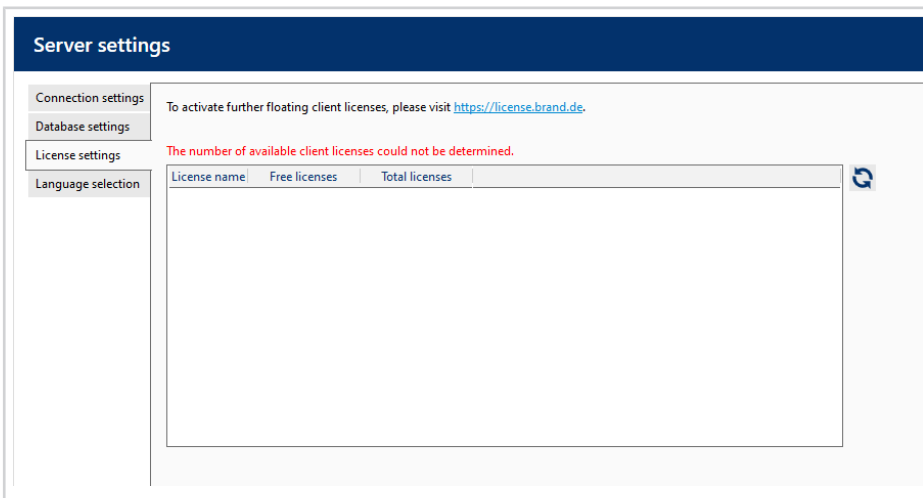
For simple networks (or if there is not already an MS SQL Server) leave the settings on SQLite. The path to the EASYCAL-database should not be changed. Moving to a "network drive" is not supported; paths such as [\\directory\network](#) are not supported.

If you want to use a different folder on a **local drive**, make sure that the folder was created by the software. Otherwise, SQLite database `easyCalServer.db` is not created.

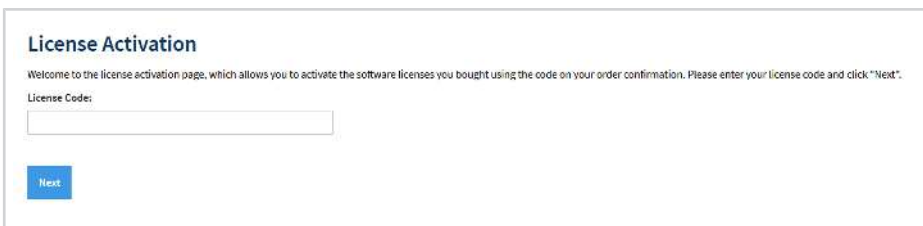
## 6 Add client licenses.

- a. Go to tab License settings.

- b. Click on the URL in the window, the default browser will open.

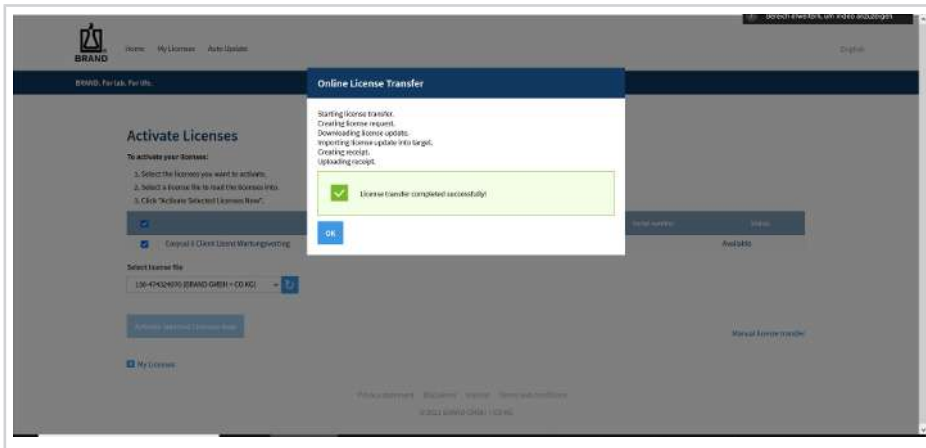
- c. Follow the license activation dialog analogously to the license activation of the server described above. Client licenses can be purchased via the BRAND Webshop (<https://shop.brand.de>).
- d. Select the license container of the server. Only one container should be available.

- e. Click on Activate Selected Licenses Now .

↪ The license is now activated.

Add client licenses.

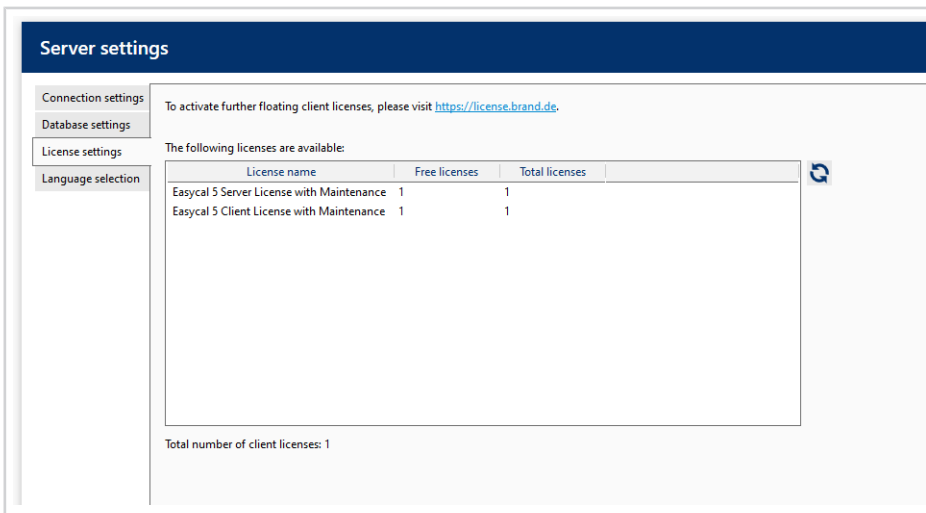
f. Confirm with OK.

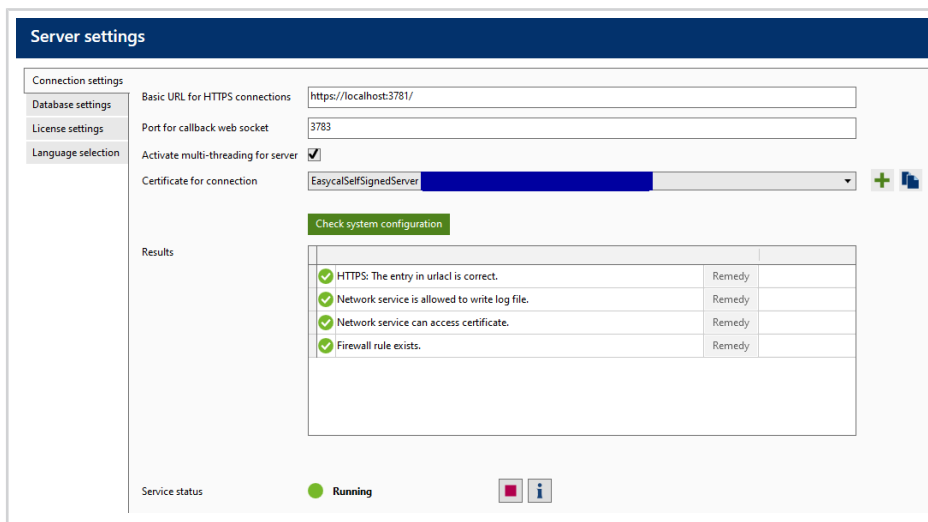
g. After completing the activation in the web browser, switch back to the EASYCAL™ 5 server software.

h. In the License settings tab click the button .

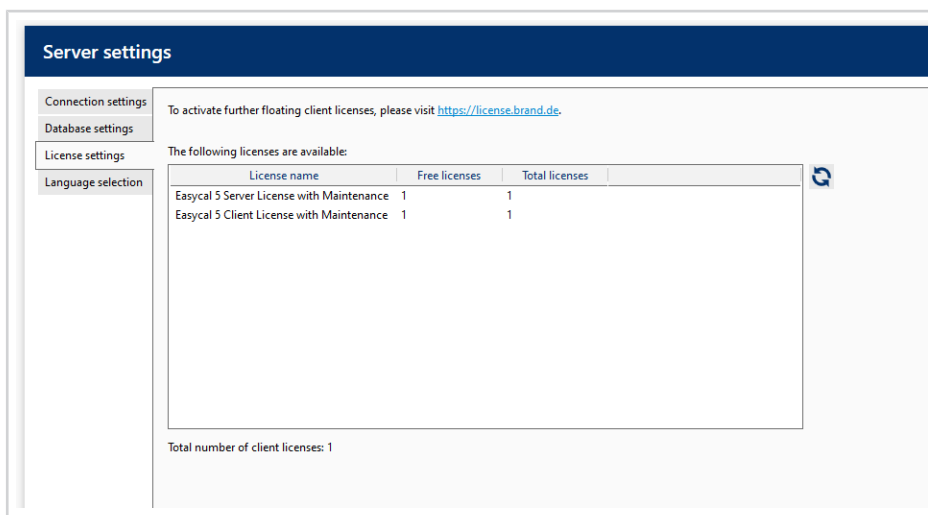
i. In this way, any number of clients can be linked to the server.

j. Finally, switch to the Connections settings tab and configure the server by clicking the button .

➤ This completes the preparations on the server.

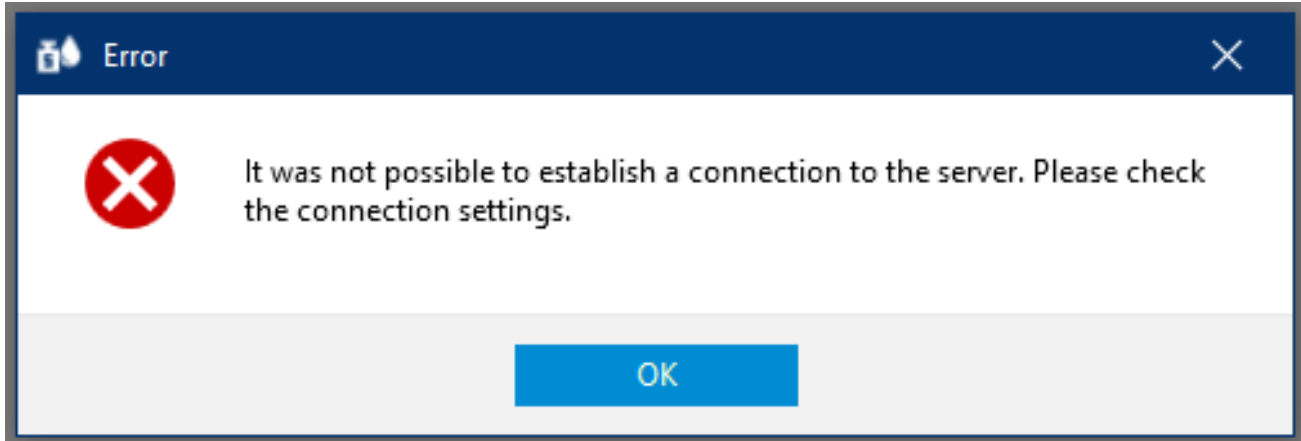
- k. From now on, the EASYCAL™ 5 client software can be installed on the client PCs in the network. As many client PCs with EASYCAL™ 5 can be operated in parallel as there are unused client licenses available on the EASYCAL™ 5 server. A current overview can be seen in the License settings tab.

## 7 First Client Start

a. Start the client with local admin rights.

→ After the start an error message appears, because the EASYCAL Client Software does not know the address to the EASYCAL server yet.

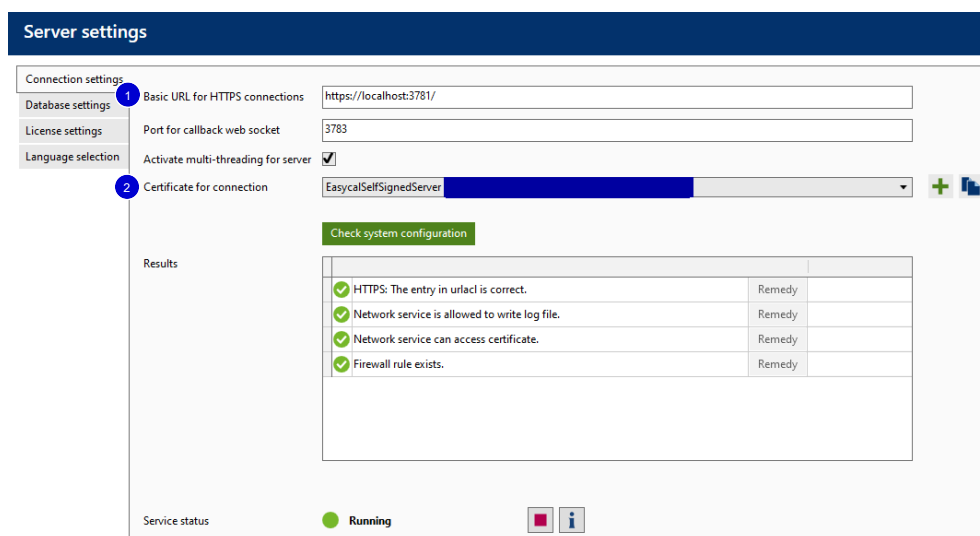
→ If the EASYCAL server is already configured and the Product activation required window appears when the client is started, the EASYCAL client cannot connect to the license server over the network. Remedy, see Configure the license server on the client, p. 41

b. Confirm Error message by clicking OK.

→ The window for entering the EASYCAL server address appears.

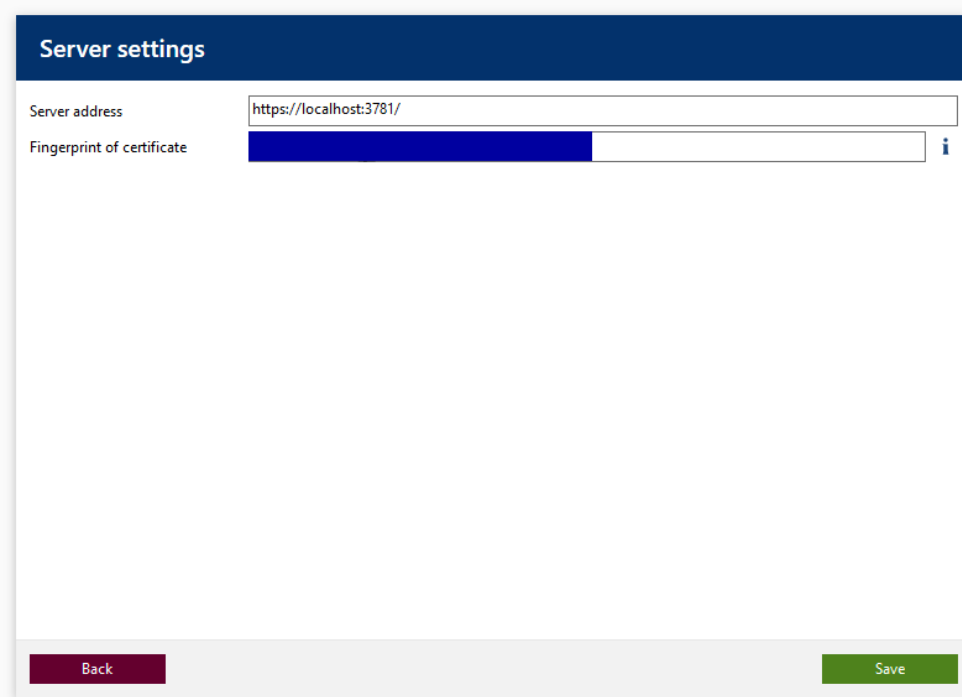
c. Copy the Basic URL and the Certificate for connection from the server into a text file. Then insert the entries as follows in the Client Software > Settings > Server settings.

Basic URL for HTTPS connections	insert in	Server address
Certificate for connection	insert in	Fingerprint of certificate

1 Basic URL/Server address

2 Certificate for connection/finger print

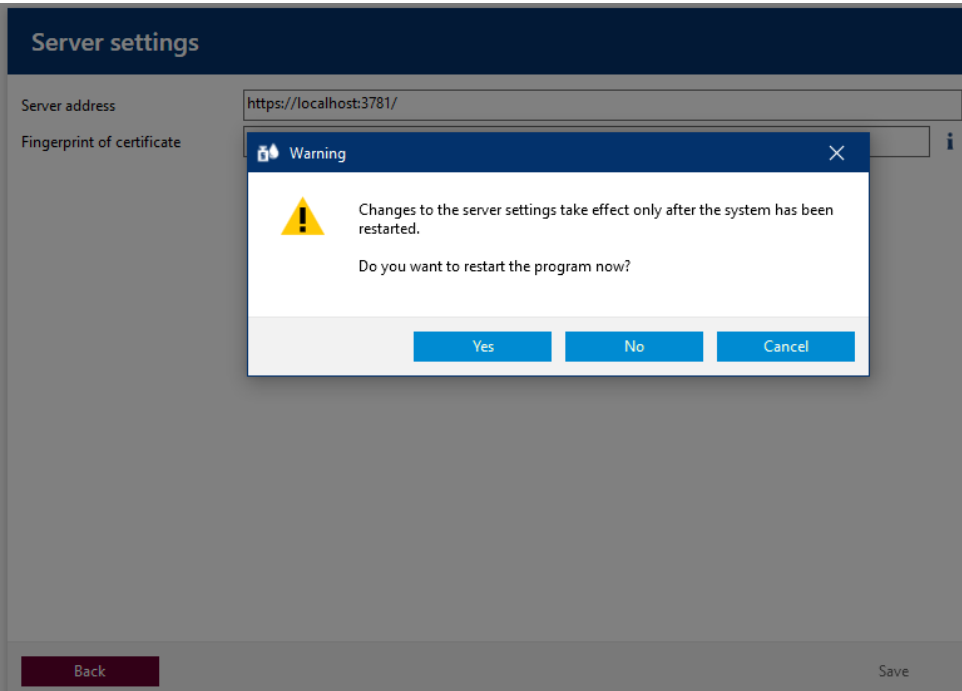
**Server settings**

Server address

Fingerprint of certificate

Back Save

d. Click on Save in the EASYCAL client software.

**Server settings**

Server address

Fingerprint of certificate

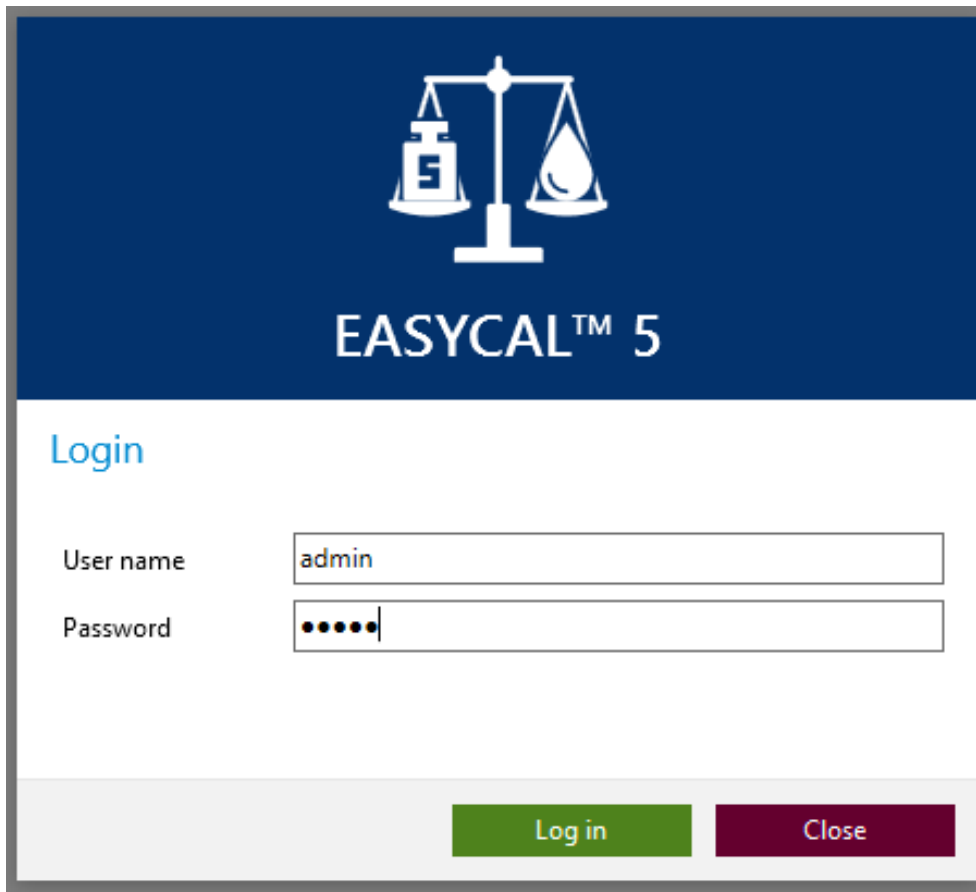
Warning

⚠ Changes to the server settings take effect only after the system has been restarted.  
Do you want to restart the program now?

Yes No Cancel

Back Save

e. Restart EASYCAL client.

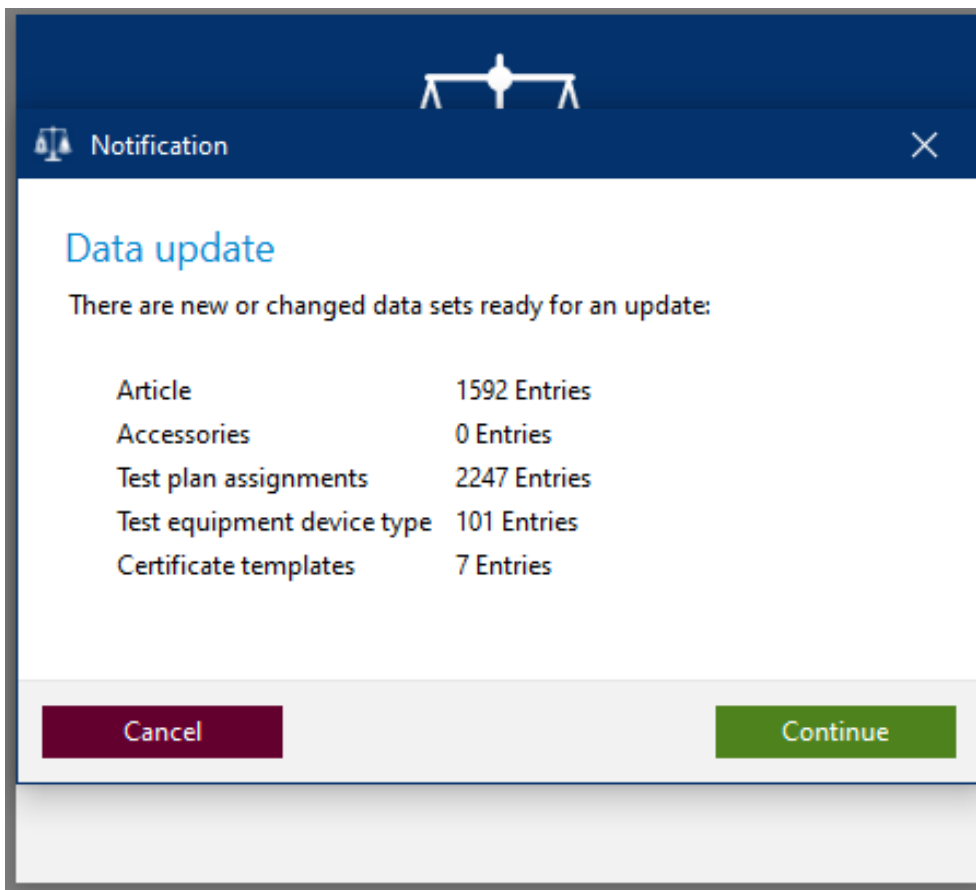
**EASYCAL™ 5**

Login

User name

Password

- f. Login with standard user (user: admin | password: start) in the EASYCAL client. If necessary, an update is installed, e.g. during the initial installation.

g. After this is finished, further EASYCAL clients can be started - if free licenses are available.

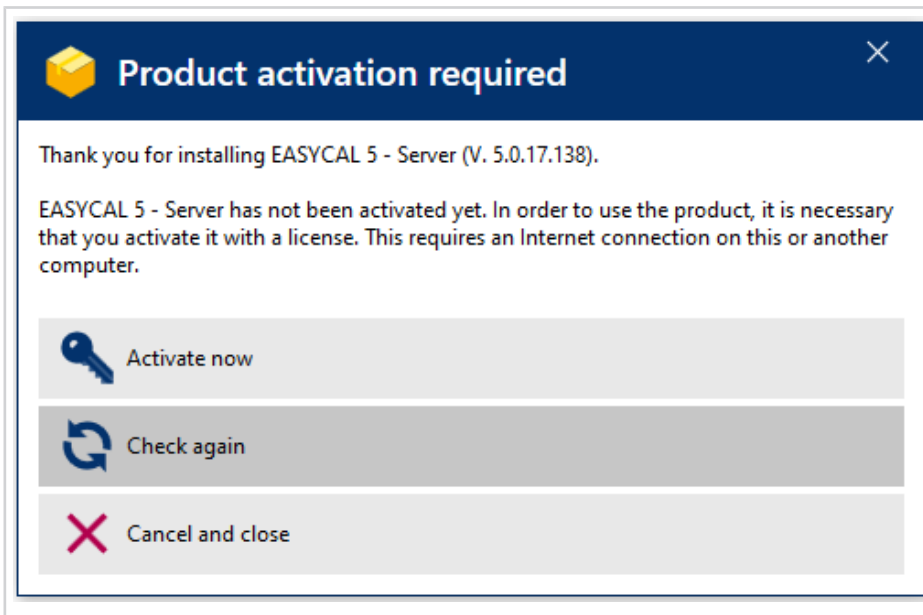
### NOTICE

#### Change default user

After completing the configuration work, change the default user (user: admin | password: start ) in order to prevent unauthorized access to the system.

#### EASYCAL client does not start | Product activation required appears

The EASYCAL client does not start even though the connection to the EASYCAL server has been configured. The Product activation required window appears because EASYCAL cannot reach the license server (network structure, security settings). To ensure that EASYCAL can still be operated, the license server can be accessed via the CodeMeter Control Center , see Configure the license server on the client, p. 41.

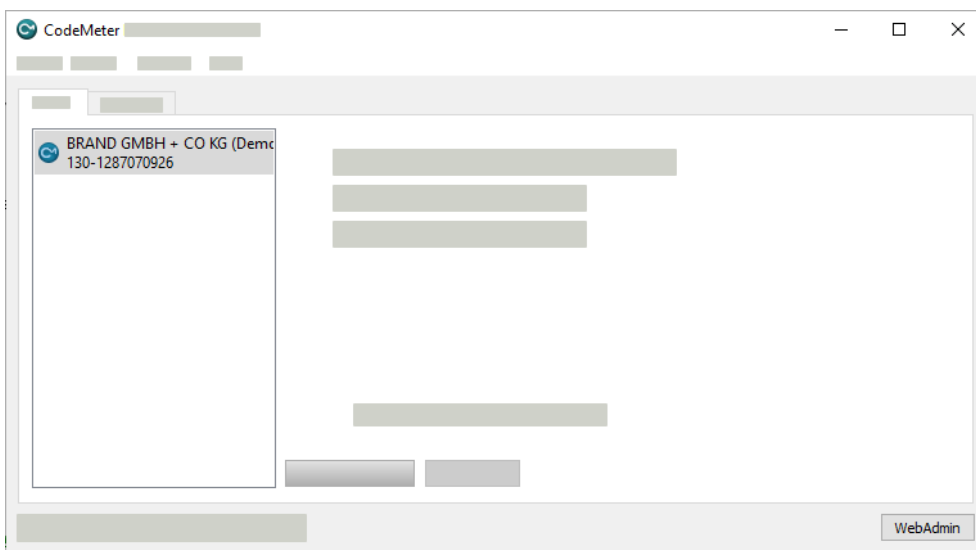
## 8 Configure the license server on the client

The license server can also be configured on the client via the CodeMeter Control Center if other connections to the license server over the network fail or are not possible.

Prerequisite:

- ✓ The IP address of the EASYCAL server must be known. Example: 192.168.3.169

- EASYCAL™ 5 Close.
- Open the CodeMeter Control Center and use the WebAdmin button to open the Web Interface.

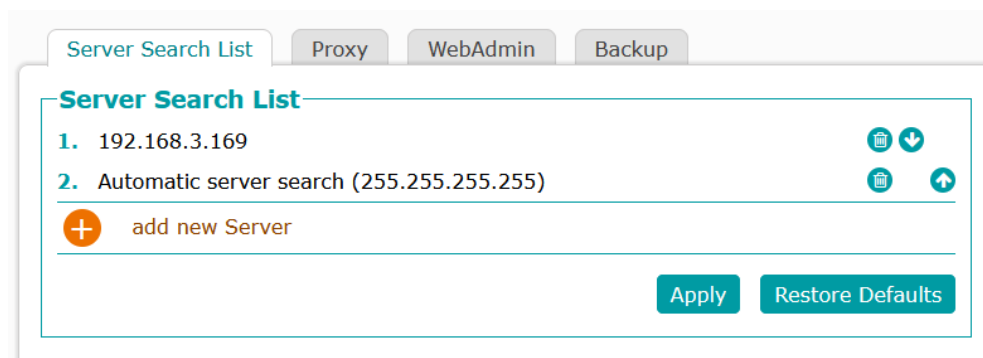
- From there, open Settings > Base > Server Search List in the navigation bar.

- Click add new Server .
- Enter the IP address 192.xxx.xxx.xxx of the EASYCAL server in the field and accept the entry. Example: 192.168.3.169

Configure the license server on the client

f. Start EASYCAL™ 5.